

Technical Note Buffer dispensing from jerrycans

Product Description

For convenient use of small volumes of buffers from the bulk supply in jerrycans we provide a set of reusable dispensing taps. The taps allows filling of pipetting reservoirs or smaller containers for intermediate use of the buffers depending on the customer's workflow requirements.

Important user notes:

- Use only 1 dispensing tap per buffer type. Contamination from 1 both buffer to another may result in quality issues (e.g. salt contamination in the Elution Buffer)
- After use, wipe the tap with a fiber-free cloth to prevent spillage

Use of the dispenser taps:

1) Remove the closure lid of the jerrycan



2) Screw the dispensing tab onto the jerrycan



3) Place the jerrycan on its side and open the dispensing tab to transfer the buffer solution to a container of your preference



Table 1: Reagents and equipment

Product	Art. No.	Contents
MagSi-NA Pathogens	MDKT0021025K	25000 preps
Dispenser Tap set for use with jerry cans	MDPL00270005	Pack of 5 pieces

magtivio B.V.
Daelderweg 9
6361 HK Nuth (The Netherlands)

Phone: +31-(0)45-2084810
Fax: +31-(0)45-2084817
Internet: www.magtivio.com