

Technical Note

Buffer dispensing from jerrycans

Product Description

For convenient use of small volumes of buffers from the bulk supply in jerrycans we provide a set of reusable dispensing taps. The taps allows filling of pipetting reservoirs or smaller containers for intermediate use of the buffers depending on the customer's workflow requirements.

Important user notes

- Use only 1 dispensing tap per buffer type. Contamination from 1 both buffer to another may result in quality issues (e.g. salt contamination in the Elution Buffer)
- After use, wipe the tap with a fiber-free cloth to prevent spillage

Use of the dispenser taps

- 1) Remove the closure lid of the jerrycan



- 2) Screw the dispensing tab onto the jerrycan



- 3) Place the jerrycan on its side and open the dispensing tab to transfer the buffer solution to a container of your preference



Table 1. Reagents and equipment

Product	Art. No.	Contents
MagSi-NA Pathogens	MDKT0021025K	25000 preps
Dispenser Tap set for use with jerrycans	MDPL00270005	Pack of 5 pieces